


QUICK GUIDE

Monkey NGAL ELISA Kit

KIT 045

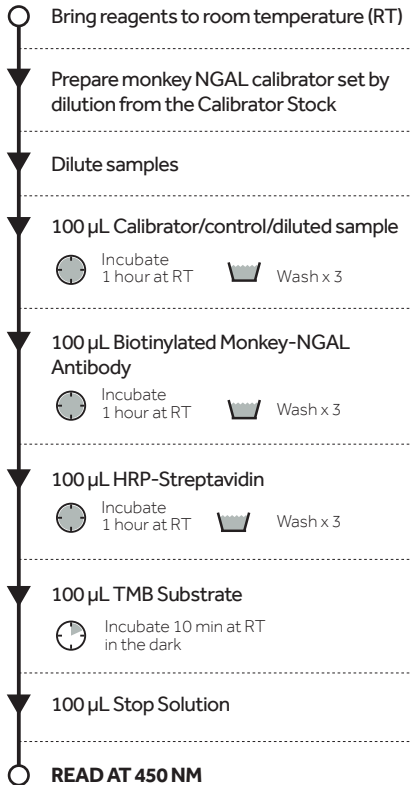


bioporto


DIAGNOSTICS

Phone (+45) 4529 0000 | info@bioporto.com | www.bioporto.com

Calibrator dilution guide*

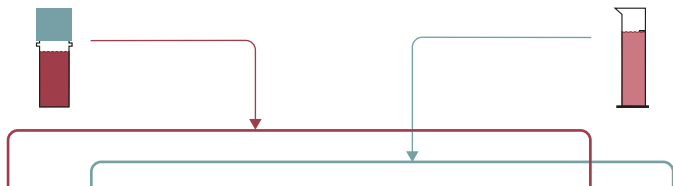

| Dilution Factor | 1 | 2 | 4 | 10 | 20 | 40 | 100 | - |
|-----------------|--------|--------|--------|--------|--------|--------|--------|--------|
| Monkey NGAL | 200 µL | 200 µL | 100 µL | 40 µL | 20 µL | 10 µL | 10 µL | - |
| Sample Diluent | - | 200 µL | 300 µL | 360 µL | 380 µL | 390 µL | 990 µL | 400 µL |

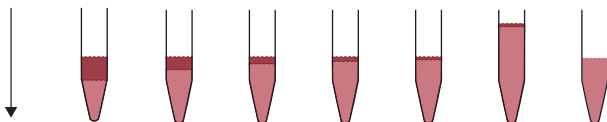

| Calibrator # | 8 | 7 | 6 | 5 | 4 | 3 | 2 | 1 |
|--------------------|------------|-----------|-----------|-----------|----------|----------|----------|---------|
| NGAL Concentration | 1000 pg/mL | 500 pg/mL | 250 pg/mL | 100 pg/mL | 50 pg/mL | 25 pg/mL | 10 pg/mL | 0 pg/mL |

| | | | | | | | | |
|-------------------|---------------------|---------------------|---------------------|---------------------|---------------------|---------------------|---------------------|---------------------|
| Transfer to wells | 100 µL/ well x 2 | 100 µL/ well x 2 | 100 µL/ well x 2 | 100 µL/ well x 2 | 100 µL/ well x 2 | 100 µL/ well x 2 | 100 µL/ well x 2 | 100 µL/ well x 2 |
|-------------------|---------------------|---------------------|---------------------|---------------------|---------------------|---------------------|---------------------|---------------------|

*Calibrators are to be freshly diluted immediately prior to use.

