


QUICK GUIDE

Monkey NGAL ELISA Kit

REF KIT 045

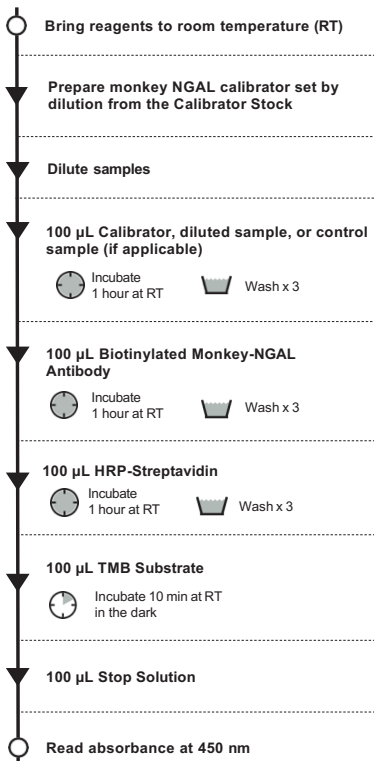

bioporto


DIAGNOSTICS

Phone (+45) 4529 0000 | info@bioporto.com | www.bioporto.com


Calibrator dilution guide*

Monkey NGAL Calibrator Stock

1x Sample Diluent

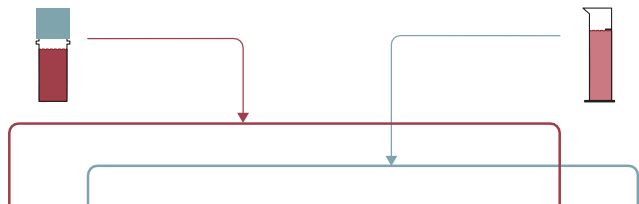

Dilution Factor	1	2	4	10	20	40	100	-
Monkey NGAL	200 μ L	200 μ L	100 μ L	40 μ L	20 μ L	10 μ L	10 μ L	-
Sample Diluent	-	200 μ L	300 μ L	360 μ L	380 μ L	390 μ L	990 μ L	400 μ L

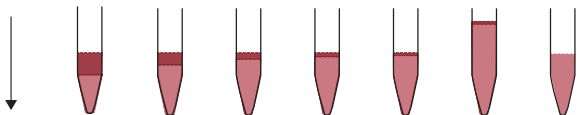

Calibrator #	8	7	6	5	4	3	2	1
NGAL Concentration	1000 pg/mL	500 pg/mL	250 pg/mL	100 pg/mL	50 pg/mL	25 pg/mL	10 pg/mL	0 pg/mL

Transfer to wells	100 μ L/ well x 2	100 μ L/ well x 2	100 μ L/ well x 2	100 μ L/ well x 2	100 μ L/ well x 2	100 μ L/ well x 2	100 μ L/ well x 2	100 μ L/ well x 2
-------------------	--------------------------	--------------------------	--------------------------	--------------------------	--------------------------	--------------------------	--------------------------	--------------------------

*Calibrators are to be freshly diluted immediately prior to use.

